Inspired by **temperature**



Connectivity

Data communication solutions for temperature control devices



Connectivity

Introduction



Table of contents

| Overview matrix | 4 - 5 |
|--|--------|
| Device groups, interfaces, protocols | |
| Available parameters | 6 - 7 |
| Only valid for devices with Pilot ONE, depending on the E-grade® | |
| Interfaces & protocols | 8 - 21 |
| Ethernet RJ45 | |
| RS232 | |
| Mini-USB data communication | |
| USB-RS232 Converter | |
| ComG@te (LEVEL / POKO / AIF / ECS / RS232) | |
| Profibusgateway 2nd/3rd Edition | |
| Profinet-Gateway | |
| EtherCAT-Gateway | |
| Software Data communication | |
| SpyControl | |
| Pilot Remote Software / PyCommand | |
| Connection / Accessories | |
| Remote operation standard variant | |
| Remote operation example variant | |

Overview matrix

Device groups, interfaces, protocols

| Controller | Device series | Interface included | Interface optional | Protocol | |
|------------|-----------------------------|--------------------|--|--------------------|--|
| | | RS232 9 pin Sub-D | - | Huber: PP, LAI | |
| KISS | Immersion circulators | Mini-USB | - | Huber: PP, LAI | |
| | | - | Pt100 sensor connection optional (only at the factory) | none | |
| | | RS232 9 pin Sub-D | - | Huber: PP, LAI | |
| | | Mini-USB | - | Huber: PP, LAI | |
| OLÉ | Minichiller Unichiller | _ | 2. RS232 9 pin Sub-D Optional (only at the factory) | Huber: PP, LAI | |
| | | - | POKO / ECS Optional (only at the factory) | none | |
| | | - | Pt100 sensor connection Optional (only at the factory) | none | |
| Pilot ONE | | | - | Huber: PB, PP, LAI | |
| THOUGHE | all | RS232 15 pin Sub-D | Profibus | Profibus DP-V0 | |
| | | | Honous | Profibus DP-V1 | |
| | Ministat, CC, Unichiller | RS232 9 pin Sub-D | - | Huber: PB, PP, LAI | |
| | | | | Huber: PB, PP, LAI | |
| Pilot ONE | alle | RJ45 Ethernet | | Modbus TCP | |
| FILOUNE | ane | NJ4J LUIEITIEL | _ | OPC UA | |
| | | | | Profinet | |
| | | | | EtherCAT | |
| Pilot ONE | alle | USB | - | Huber: PB, PP, LAI | |
| Pilot ONE | alle | _ | Analogue interfaces 1 input, 3 outputs 4 mA - 20 mA / 0 V - 10 V | none | |
| Pilot ONE | alle | - | Digital interfaces POKO / ECS | none | |

| E-grade | Basic | Exclusive | Professional | Explore | DV-E-grade | OPC UA |
|---------|----------|-----------|--------------|---------|------------|--------|
| Cat.No. | Standard | #9495 | #9496 | #10495 | #9943 | #10561 |

| Additionally required hardware | | Required E-grade level |
|---|---------|------------------------|
| Designation | Cat.No. | Designation |
| none | - | none |
| none | - | none |
| Pt100 sensor connection for KISS (only measurement, no control) | #10688 | none |
| none | - | none |
| none | - | none |
| 2. RS232 interface OLÉ | #10726 | none |
| Interface module for units with OLÉ controller | #10689 | none |
| Pt100 sensor connection for OLÉ controller (only measuring, no control) | #10519 | none |
| none | - | from Basic |
| RS232 to Profibus DP-V0 | #522248 | from Basic |
| RS232 to Profibus DP-V1 | #10503 | from Basic |
| none | - | OPC UA |
| Ethernet to Profinet Gateway | #10965 | from Basic |
| Ethernet to EtherCAT Gateway | #10966 | from Basic |
| - | - | from Basic |
| Com.G@te external for Unistats | #6915 | none |
| Com.G@te internal for all other temperature control units (except CC) | #31217 | none |
| POKO / ECS Interface for Unistats as standard 2. POKO / ECS with Com.G@te external | - | none |
| Internal POKO / ECS interface for all other temperature control units | #10003 | none |
| Com.G@te internal for all other temperature control units (except CC) | #31217 | none |

Available parameters only valid for units with Pilot ONE®

| Adress (hex) | Description | Basic Standard | Exclusive #9495 | Professional #9496 | DV / OPC UA / Explore #9943 / 10561 / 10495 |
|-----------------|--|--------------------------|---------------------------|------------------------------|---|
| 00 | Setpoint temperature controller | | \diamond | <i>🗞</i> | Ø |
| 01 | Internal temperature | | | <i>🗞</i> | Ø |
| 02 | Return flow temperature | - | - | - | Ø |
| 03 | Pump pressure (absolute) | $\boldsymbol{\diamond}$ | \diamond | ∽ | Ø |
| 04 | Current output | - | - | - | Ø |
| 05 | Fault report | \diamond | \diamond | \diamond | Ø |
| 06 | Warning report | \$ | \diamond | \diamond | Ø |
| 07 | Process temperature (Lemosa) | \diamond | \diamond | <i>🗞</i> | ∽ |
| 08 | Actual value specification internal temperature | - | - | - | Ø |
| 09 | Specification process temperature | - | - | - | 8 |
| 0A | Status of the thermostat | Ø | \diamond | <i>🗞</i> | Ø |
| OB | Activate the Blow-Down slider | \diamond | \diamond | <i>🗞</i> | Ø |
| 0C | Activate Blow-Down slider heating | \diamond | \diamond | <i>🗞</i> | Ø |
| OF | Fill level | \diamond | \diamond | <i>~</i> | Ø |
| 12 | PID parameter automatic temperature controller | \diamond | Ø | <i>~</i> | Ø |
| 13 | Temperature control mode | - | \diamond | <i>∞</i> | Ø |
| 14 | Temperature control | \$ | Ø | <i>~</i> | Ø |
| 15 | Operating mode compressor | \diamond | \diamond | <i>~</i> | ∽ |
| 16 | Circulation | \$ | <i>\lambda</i> | <i>~</i> | Ø |
| 17 | Operating lock | \diamond | \diamond | <i>~</i> | ∽ |
| 18 | Mode actual value specification internal temperature | - | - | - | |
| 19 | Mode actual value specification process temperature | - | - | - | < |
| 1A | Freeze protection | <i>⊗</i> | <i>⊗</i> | <i>⊗</i> | < |
| 1B + 1C | Serial number | < | <i>⊗</i> | <i>⊗</i> | < |
| 1D | Kp of the internal controller | <i>⊗</i> | <i>⊗</i> | <i>⊗</i> | < |
| 1E | Th of the internal controller | ✓ | <i>⊗</i> | <i>⊗</i> | × |
| 1F | Tv of the internal controller | \$ | <i>⊗</i> | | |
| 20 | Kp of the jacket controller | - | | | × |
| 21 | Tn of the jacket controller | - | <i>⊗</i> | | ✓ |
| 22 | Tv of the jacket controller | - | | | × |
| 23 | Kp of the process controller Tn of the process controller | - | <i>⊗</i> | | ✓ ✓ |
| 24 | Tv of the process controller | - | ~ ⊘ | | ✓ |
| 25 | Pump speed | ~ | | | × * |
| 20 2C | Cooling water inlet temperature | ~ | ~ | - | × ⊗ |
| 2C 2D | Cooling water inter temperature | _ | _ | _ | × |
| 2D 2E | Network connection conditions | _ | - | _ | ~ ⊘ |
| 30 | Minimum setpoint | \$ | \$ | \$ | |
| 31 | Maximum setpoint | ~ ∽ | ~ ~ | ~ ~ | ✓ |
| 33 | Over-level limit | × | ~ ~ | × | |
| 34 | Sub-level limit | × × | ~ ~ | ~ ~ | ~ ⊘ |
| 35 | Setting the digital level output | × | \$ \$ | \$ \$ | Ŷ |
| 35 3A | Process temperature | _ | ~ ⊘ | < | ✓ |
| 3C | Status of the thermostat | \$ | <i>⊗</i> | <i>⊗</i> | ✓ |
| 3D | Interference variables | - | - | - | ✓ |

depending on the E-grade®

| Adress (hex) | Description | Basic Standard | Exclusive #9495 | Professional #9496 | DV / OPC UA / Explore #9943 / 10561 / 10495 |
|-----------------|--|--------------------------|---------------------------|------------------------------|---|
| 3E | Pressure in return (absolute) | \$ | \$ | <i>~</i> | \$ |
| 3F | Status Blow-Down | \diamond | \$ | <i>🕫</i> | <i>🗞</i> |
| 40 | Watchdog (Fault) | \$ | \diamond | <i>~</i> | Ø |
| 41 | Watchdog (2nd setpoint) | - | - | <i>🗞</i> | Ø |
| 42 | 2. Setpoint | - | - | <i>🗞</i> | Ø |
| 43 | PMA Mode | - | - | - | < |
| 44 | Specify PMA output | - | - | - | ∽ |
| 48 | Setpoint pump speed | \diamond | \diamond | <i>~</i> | Ø |
| 49 | Setpoint pump pressure | \$ | <i>\U</i> | <i>~</i> | Ø |
| 4A | VPC bypass operating mode | \diamond | \diamond | <i>~</i> | Ø |
| 4B | VPC bypass target position | \$ | <i>⊗</i> | <i>~</i> | Ø |
| 4C | Cooling water outlet temperature | - | - | - | Ø |
| 4D | Thermal fluid volume flow | - | - | - | Ø |
| 4E | Setpoint thermal fluid volume flow | - | - | - | ∽ |
| 4F | Setpoint Delta-T control | - | <i>⊗</i> | <i>~</i> | ∽ |
| 50 | Alarm limit Delta-T | - | \diamond | \$ | Ø |
| 51 | Upper alarm limit internal temperature | \$ | ✓ | <i>~</i> | Ø |
| 52 | Lower alarm limit internal temperature | \$ | \diamond | <i>~</i> | < |
| 53 | Upper alarm limit process temperature | ✓ | × | <i>~</i> | ✓ |
| 54 | Lower alarm limit process temperature | ✓ | <i>~</i> | | <i>⊗</i> |
| 55 | Setting of heating overtemperature protection | ✓ | × | <i>~</i> | < |
| 56 | Setting of overtemperature protection expansion vessel | \$ | <i>~</i> | | Ø |
| 58 | Start temperature control program | - | <i>⊗</i> | <i>⊗</i> | < |
| 59 | Specify ramp duration | - | <i>⊗</i> | <i>⊗</i> | < |
| 5A | Start ramp | - | <i>⊗</i> | <i>⊗</i> | < |
| 5B | Specify Blow-Down operating status | <i>⊗</i> | <i>⊗</i> | <i>⊗</i> | <i>∞</i> |
| 5C | Days until maintenance | <i>\U</i> | <i>⊗</i> | <i>⊗</i> | < |
| 5D | Days to F-Gas check | <i>⊗</i> | <i>⊗</i> | <i>⊗</i> | <i>∞</i> |
| 5E | Create service packet | Ø | <i>⊗</i> | | < |
| 5F | Change program status | - | | | <i>∞</i> |
| 62 | Pressure of the VPC bypass | <i>\</i> | <i>~</i> | <i>⊗</i> | ✓ |
| 69 | Mode actual value specification thermal fluid volume flow | - | - | - | ♦ |
| 6A 6B | Specification thermal fluid volume flow Control mode of the pump | - | - | - | ✓ ✓ |
| 6C | External POKO control | ~ | ~ | ~ | ♥ |
| 6D | POKO status | - | - | _ | ♥ |
| 6E | Current output (high bytes) | - | - | _ | ♥ |
| | Ventilation | ~ | - | ~ | ✓ |
| 6F 70 | Emptying | ♥ | | | ♥ |
| 70 | Setpoint temperature controller | ✓ | ✓ | | ✓ |
| 72 | VPC bypass position | ✓ | ~ ⊘ | | × * |
| 72 | Message | ~ ∽ | ✓ | ✓ | ✓ |
| 73 | Disturbance feedforward VPC | _ | _ | _ | ✓ |
| 75 | Actual value specification mode pump pressure | _ | _ | | ✓ |
| 76 | Actual value specification pump pressure (relative) for pump control | | _ | _ | ~ ⊘ |
| 70 | netual value specification pump pressure (relative) for pump control | - | - | _ | Ÿ |

*same meaning / function as address 0. Is required for some OPC UA applications.


Interfaces RJ45 Ethernet (RJ45)

Standard LAN connection, no driver required.


The following components are used for an Ethernet setup:

- Cable with RJ45 connector
- Commercially available switch -> when connecting several Pilot ONEs

TCP-IP Setup Modbus TCP Setup OPC UA Setup

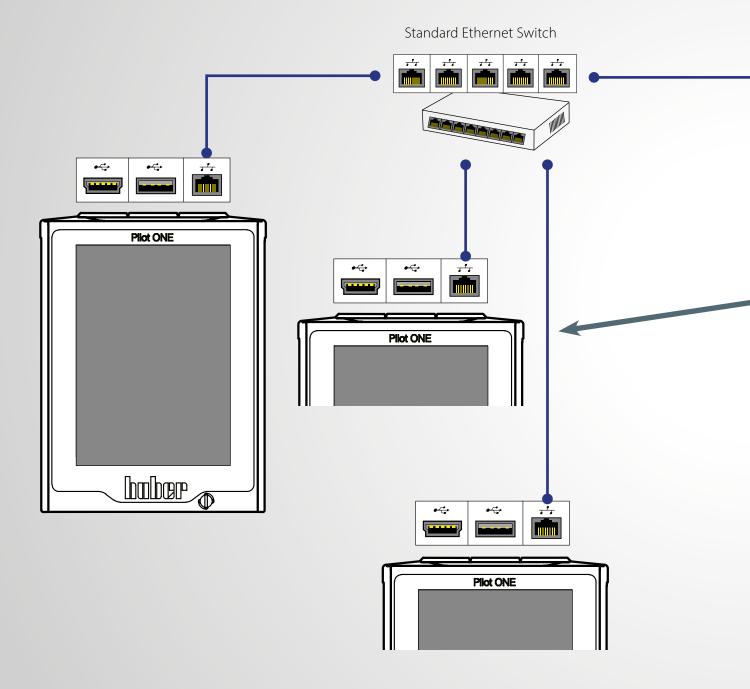

only for units with $\mathsf{Pilot}\ \mathsf{ONE}^{\texttt{B}}$

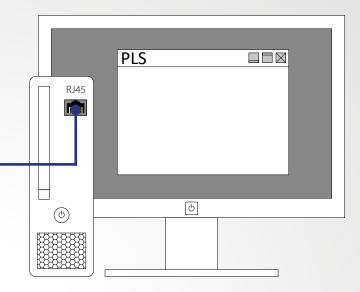

Computer or PLC with Ethernet interface

| Bus system: Cable type: | Fast Ethernet 10/100 Standard Ethernet100 Base-TX CAT5 10/100 RJ45 |
|----------------------------|--|
| Communication via: | TCP (Transmission Control Protocol) ➤ Port 8101 Modbus TCP ➤ Port 502 OPC UA ➤ Port 4840 |

| Protocol | Additionally required hardware | Additional E-grade | Cat.No. |
|-------------|--------------------------------|-----------------------|--------------|
| PB, PP, LAI | none | command dependent | see page 6/7 |
| Modbus TCP | none | command dependent | see page 6/7 |
| OPC UA | none | OPC UA E-grade | #10561 |


9

Interfaces **RS232**

Pilot ONE®-Controller


Known standard interface Point-to-point connection

Various Sub-D connectors: 15 pin on the service socket or 9 pin on the ComG@te

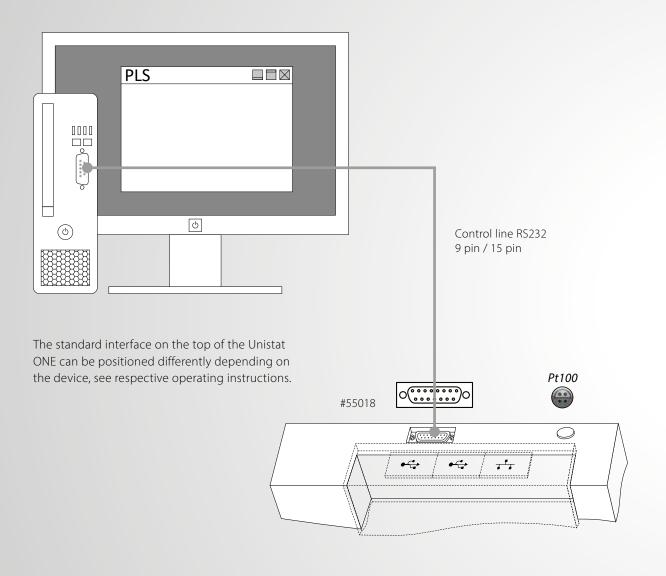


| Protocol | Controller | Socket type | Additional hardware | Additional E-grade |
|--------------------|------------|----------------|------------------------|-----------------------|
| Huber: PP, LAI | KISS, OLÉ | 9 polig Sub-D* | none | none |
| Huber: PB, PP, LAI | Pilot ONE | 15 polig Sub-D | none | command dependent |
| Huber: PB, PP, LAI | Pilot ONE | 9 polig Sub-D* | none | command dependent |
| *unit dependent | | | | |

KISS®-/OLÉ-Controller

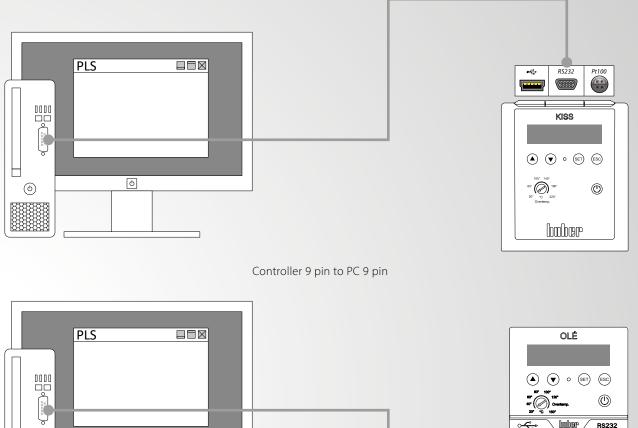

| () | ð | |
|----|---|--|
| | | |
| | | |

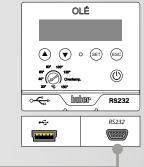


| Control cable | Length | Cat.No. |
|---|--------|----------|
| | 3 m | 55018 |
| Serial data transfer from device 15 pin Sub-D connector to PC / Notebook 9 pin Sub-D socket | 5 m | 55018-5 |
| to rey notebooks pinsus a societ | 10 m | 55018-10 |
| Serial data transmission from device / | 3 m | 6146 |
| Com.G@te 9 pin Sub-D connector to PC / Notebook 9 pin Sub-D socket | 5 m | 6146-5 |
| | 10 m | 6146-10 |



Interfaces Mini-USB data communication

With a direct USB connection, a driver must be installed for the Pilot ONE or for KISS/OLÉ that installs a virtual COM port on a Windows PC. This virtual COM port behaves like an RS232 port and is selected accordingly in communication software (e.g. SpyControl) also as an RS232 port (e.g. COM 5).

You can find the drivers on our website in the Service & Download --> Software area.

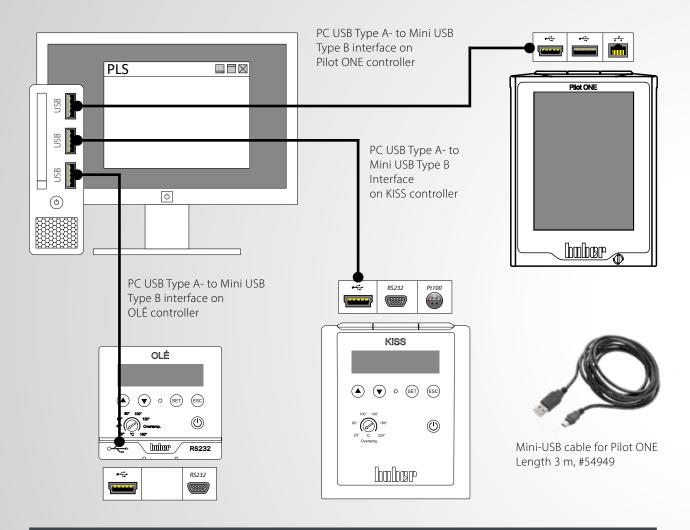

| Protocol | Controller | Socket type | Additional hardware | Additional E-grade |
|--------------------|------------|-------------|------------------------|-----------------------|
| Huber: PP, LAI | KISS, OLÉ | Mini-USB | none | none |
| Huber: PB, PP, LAI | Pilot ONE | Mini-USB | none | command dependent |


| Control cable | Length | Cat.No. |
|---|--------|---------|
| Serial data transfer from controller Pilot ONE, KISS and OLÉ to PC / notebook, USB Type A | 3 m | 54949 |

USB-RS232 Converter

When using a USB-RS232 converter, a virtual COM port is also installed on Windows PCs. This virtual COM port can also be used like a normal hardware RS232 port. Drivers are available from corresponding converter manufacturer.



Important: The converter must not be connected directly onto Huber devices - neither with Pilot ONE nor with KISS/OLÉ. A null modem cable (crossed cable), e.g. # 6146 must always be used between converter and Huber device.

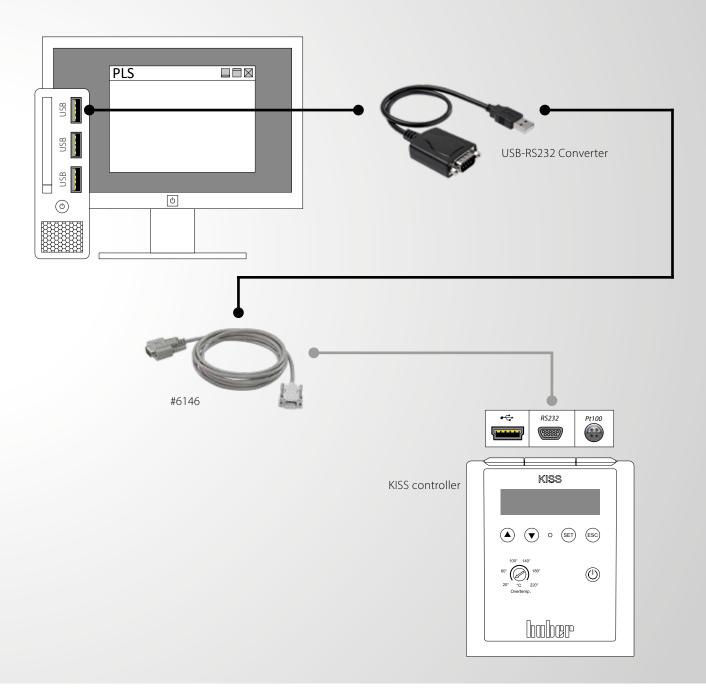


Interfaces Com.G@te

The Com.G@te provides connections according to the NAMUR standard. It expands devices with Pilot ONE with additional interfaces and connection options. The following interfaces are integrated: RS232 (bidirectional), ECS external control signal (programmable), POKO potential-free contact (programmable), AIF analogue interface 0/4-20 mA or 0-10 V (bidirectional).



Com.G@te external interfaces:

- LEVEL
- POKO
- AIF
- ECS
- RS232

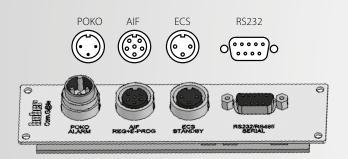

Com.G@te internal interfaces:

- POKO
- AIF
- ECS
- RS232

POKO, AIF, ECS, RS232 can be built into: Ministat, Petite Fleur, Grande Fleur, Unichiller, CCs ab CC-4xx

| Interface module | Description | Cat.No. |
|------------------|-------------|---------|
| Com.G@te intern | | 31217 |
| Com.G@te extern | | 6915 |

| Length | Cat.No. |
|--------|-------------------|
| 3 m | 9490 |
| 3 m | 9491 |
| 3 m | 9492 |
| 3 m | 9353 |
| | 3 m 3 m 3 m |

POKO connector

(Potential-free contact) alarm

Signal contact for external monitoring.

Note the functional options that POKO offers you in the "Interfaces" category. The switching behaviour can be set in the Pilot ONE, e.g. "Fault". The potential-free contact (POKO) signals the status of the temperature control unit via the contact position. A closed normally open contact means operational readiness. In the event of a fault or error, the normally open contact is opened (this applies to the closer contact between pin 1 and pin 2).

ECS

(External Control Signal) Stand-by

Control signal for activating a function. The functionality of the ECS is determined by the "Interfaces" category, e.g. the release signal ECS (external control signal) can be used to start/stop the temperature control.

The control takes place via a potential-free contact. ECS becomes electronically active when E1 and E2 are connected by an external potential-free contact. Contact specification: min. 0.1 A/24 V DC.

AIF Reg-E-Prog

The analogue interface has a programmable input channel and 3 output channels 4 - 20 mA.

The analogue interface of the Com.G@tes is programmed in the "Interfaces" category.

2 output channels are fixed:

- Output 1 Setpoint
- Output 2 Supply line temperature
- Output 3 Freely configurable

RS232

This socket offers a connection via an RS232 interface, a PLC or a distributed control system (DCS) so that remote control of the controller electronics can be connected.

When using the interface, the specifications of the generally applicable standards must be observed.

LEVEL

This connection is only available with an external Com.G@te and offers the possibility of connecting an external float switch (#6152), which is positioned in the inspection glass, for level monitoring of an externally closed application. The control takes place via a potential-free contact.

 (ι) The interface is specified as a digital input. Do not apply voltage or current!



Interfaces **Profibusgateway**



Only for units with Pilot ONE


Two versions are available:

Profibusgateway DP-V0:

Preconfigured parameter list for cyclic Profibus communication. Profibus master must support Profibus DP-V0.

Profibusgateway DP-V1:

Configurable parameters via acyclic Profibus communication. Profibus master must support Profibus DP-V1.

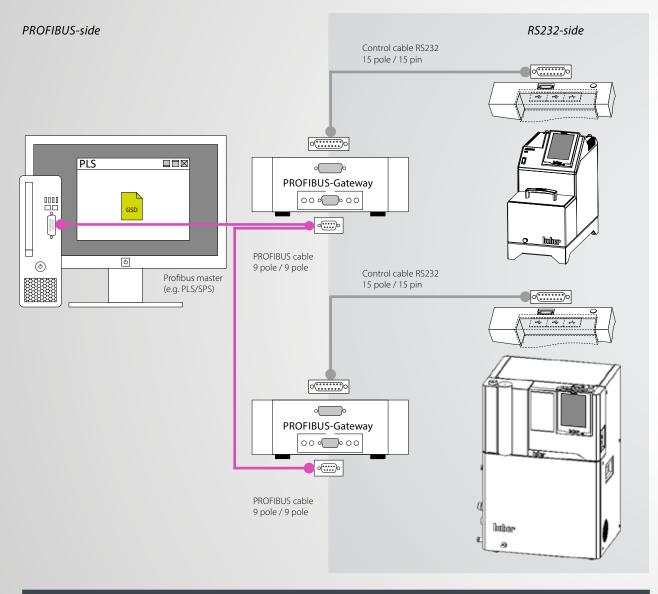


| Gateway | Description / Requirement | Cat.No. |
|------------------------|--|---------|
| Profibus Gateway DP-V0 | Parameter fix (Profibus DP-V0) | 522248 |
| Profibus Gateway DP-V1 | Parameter freely configurable (Profibus DP-V1) | 10503 |





PROFINET (Process Field Network) is the open Industrial Ethernet standard of the PROFIBUS user organisation e. V. (PNO) based on Ethernet-TCP/IP and complements PROFIBUS technology for applications that require fast data communication via Ethernet networks in combination with industrial IT functions. PROFINET can be used to implement solutions for manufacturing technology, process automation, building automation and the entire spectrum of drive technology. With the PROFINET-Gateway, Huber temperature control units can be integrated into PROFINET networks easily, flexibly and close to the process. The PROFINET-Gateway is integrated into the project planning software with the aid of the associated GSDML file.

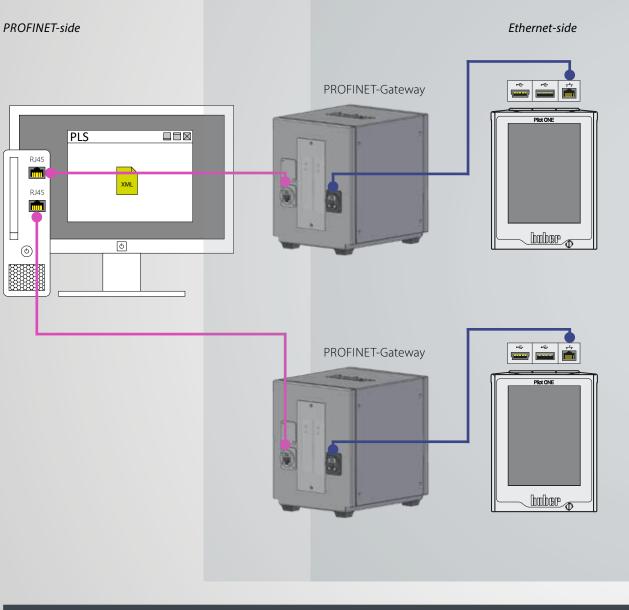


| Model | Cat.No. |
|------------------|---------|
| PROFINET-Gateway | 10965 |



Interfaces EtherCAT-Gateway EtherCAT-Gateway

EtherCAT, short for Ethernet for Control Automation Technology, is an open Ethernet-based fieldbus technology that is standardised in international standards. EtherCAT is a very fast Industrial Ethernet system that is also suitable for use in timecritical applications. With the EtherCAT gateway, Huber temperature control units can be integrated into EtherCAT networks easily, flexibly and close to the process. The EtherCAT gateway is integrated into the configuration software with the aid of the associated GSDML file.

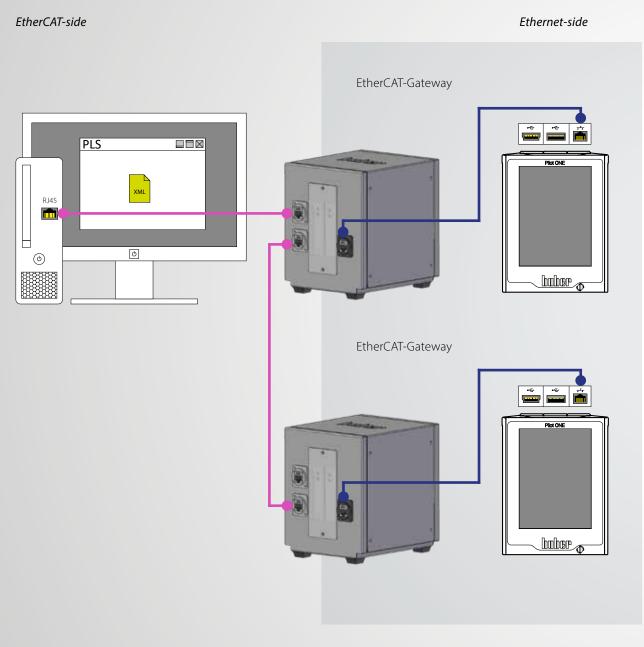

| Model | Cat.No. |
|------------------|---------|
| EtherCAT-Gateway | 10966 |



Software Data communication

<page-header><text><figure><figure>

SpyControl

SpyControl is a software solution for Windows PCs for device control as well as for the visualisation and documentation of process-relevant data. Communication with the temperature control unit takes place via RS232, USB or TCP/ IP. Recorded data are displayed over time, with the axes of the diagram being freely scalable. A zoom function simplifies the graphic evaluation of individual time segments.

The free version allows the control of a maximum of one device. The paid full version (license key required) can communicate with up to 10 devices at the same time and offers additional functions. A setpoint specification, a start/ stop function and a switchover between internal and process temperature control are possible for each channel. The

temperature values and other process data can be recorded and saved as a CSV file. Using a programmer function with a user-friendly graphic editor, an automatically running temperature profile can be specified (for max. 1 channel).

SpyControl is available as a downloaded from the Huber website. The free version is included in the download (max. 1 device) as well as a 30-day trial version of the paid full version (max. 10 devices). If you want to continue using the full version after the test period has expired, you will need to purchase a license key.

Description

SpyControl full version license key (10 channels)

66108

Cat.No.



Pilot Remote Software

The Pilot Remote software enables the complete remote control of Huber temperature control units with Pilot ONE via any PC/laptop with Microsoft Windows.

The familiar Pilot ONE user interface is displayed on the PC, i.e. identical operation on PC and temperature control unit. Communication takes place via an Ethernet network connection with secure authentication and encryption.

The software is executable under Microsoft Windows 7/8/10 (32/64 bit). To operate the software, the E-grade Remote GUI must be activated on the connected Pilot ONE units. A 30-day evaluation version can be activated directly on the Pilot ONE free of charge.

| Description | Cat.No. |
|-----------------------|---------|
| Pilot Remote Software | free |
| E-grade Remote GUI | 520450 |



PyCommand

This Python programming example makes it easier to familiarise yourself with the subject of interface communication (Ethernet, RS232, USB) with Huber temperature control devices. The components Python 2.7.4 and pySerial 2.6 from the runtime package are required.

PyCommand is a graphic program as an aid for Huber command sets. With the software, individual commands can be sent to Huber devices and commands can be tried out without having to program. The program can be used with the interfaces RS232, USB and Ethernet.


With PySoftcheck, a free Python programming example is also available as a download, which facilitates familiarization with the topic of interface communication (Ethernet, RS232, USB) with Huber temperature control units. The components Python 2.7.4 and pySerial 2.6 from the Runtime Package are required.

➡ Free download on www.huber-online.com


Connection / Accessories **Remote operation**

Standard variant

Modules that are connected as accessories to devices with Pilot ONE communicate via an internal CAN bus. For remote operation, please note that the RS232 interface can only be used in conjunction with a special cable. The interfaces on the Pilot ONE (USB / Ethernet) can be used without restrictions.

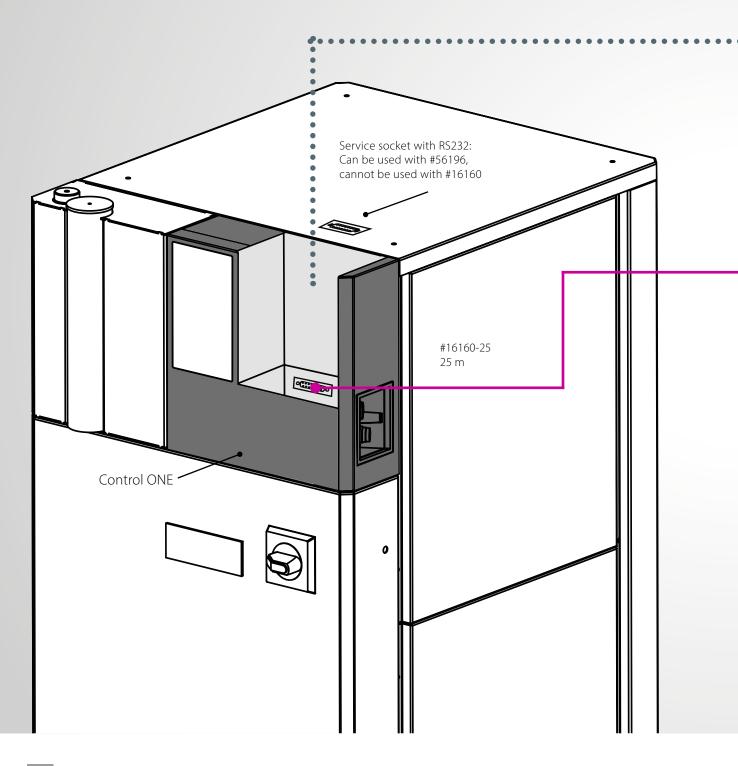

via internal CAN bus - only for units with Pilot ONE®

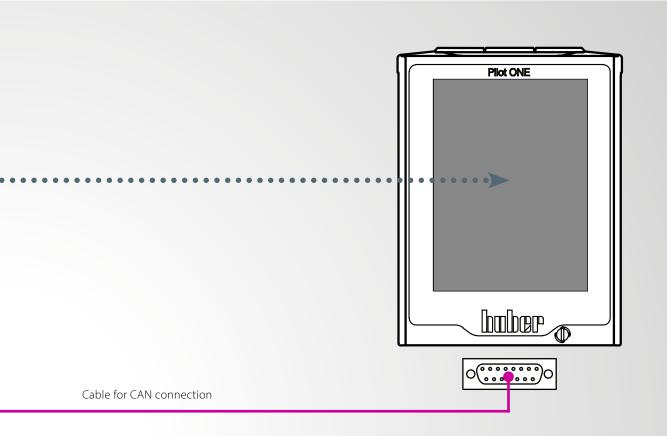

| ltem | Length | Cat.No. |
|---|---|---------|
| Extension cable for controller Pilot ONE for using the controller as remote control | 3 m (Standard) 5 m (Optional) 10 m (Optional) 15 m (Optional) 20 m (Optional) 25 m (Optional) 30 m (Optional) | 16160 |
| RS232 only possible with | 0,5 m (Optional 1 m (Optional) 3 m (Standard) 5 m (Optional) 6 m (Optional) 10 m (Optional) | 56196 |

★ Note:


If only the Pilot ONE is remote, no CAN switch is necessary. However, this can be retrofitted at any time if necessary.


25

Connection / Accessories **Remote operation**


Example variant

The Pilot ONE is used in remote operation and expanded with a Com.G@te and a VPC bypass for pressure control. A CAN switch is required for this.

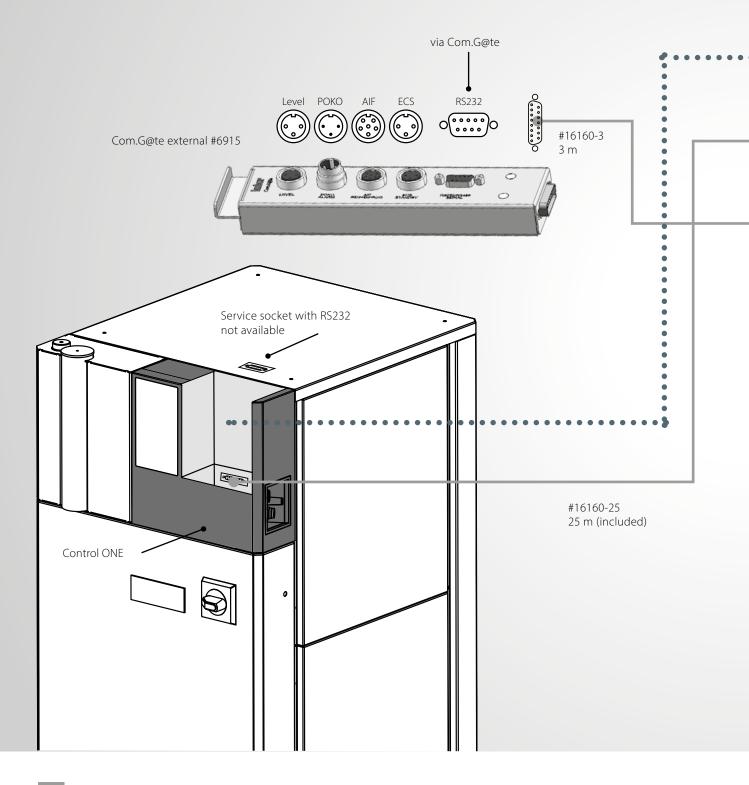

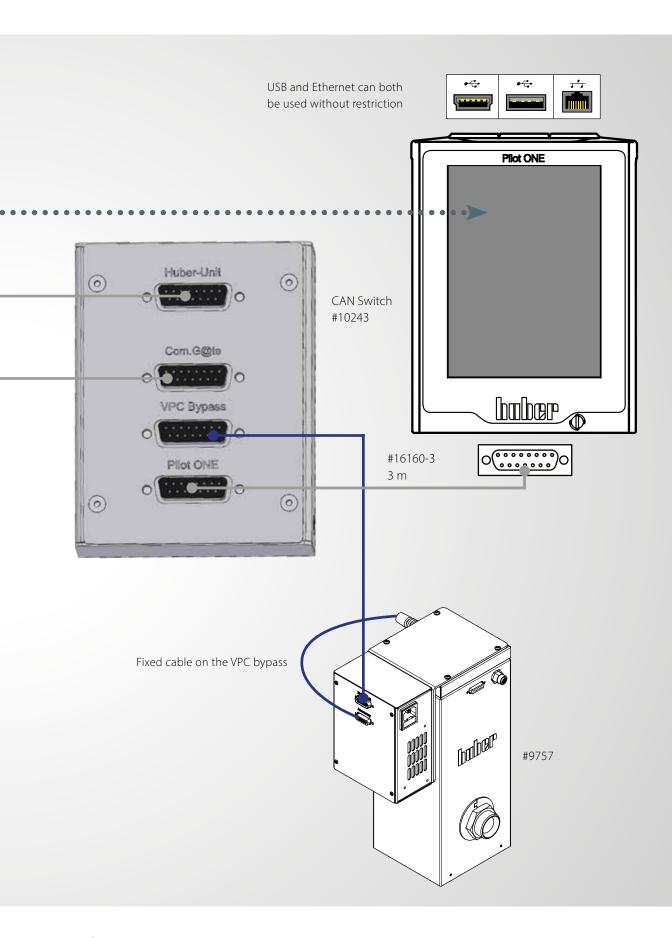

Inspired by **temperature** designed for you

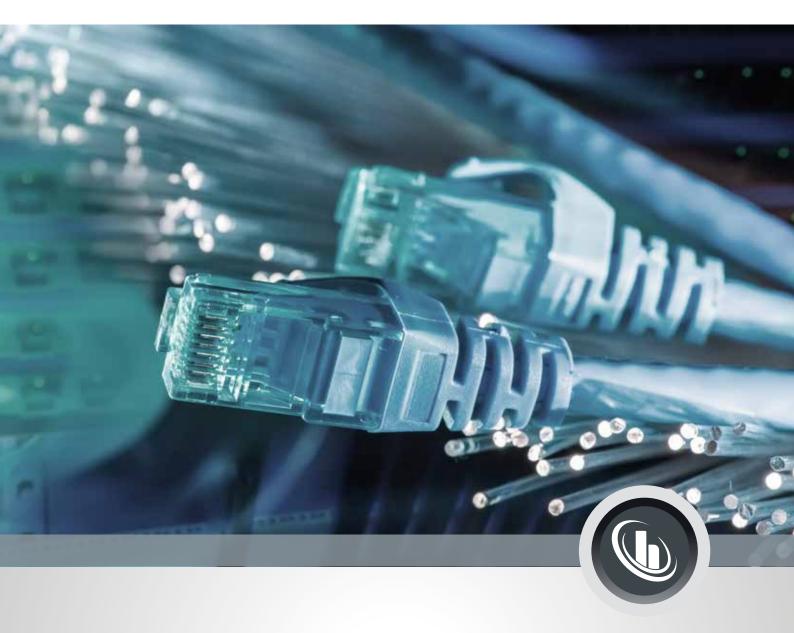

Peter Huber Kältemaschinenbau SE

Werner-von-Siemens-Str. 1 · 77656 Offenburg / Germany Phone +49 (0)781 9603-0 · Fax +49 (0)781 57211 info@huber-online.com · www.huber-online.com

